

Product number: U8476
 Product name: DCAM-API
 Product version: December 2008
 Description:

For Macintosh OSX

FireWire

DCAM Version		
DCAM Module	3.00.352.3032 2.1.3.3032	(for ORCA and C8484/C9440) (for ORCA II)
DRIVER	3.00.361.3032 2.1.3.3032	(for ORCA and C8484/C9440) (for ORCA II)
Supported Camera		
C10600-10B	ORCA-R2	
C4742-80-12AG	ORCA-AG	
C4742-95-12ERG	ORCA-ERG	
C4742-95-12NRG	ORCA-NRG	
C4742-95-12G04 / 96-12G04	ORCA-285G	(*1)
C4742-98-ERG(2)	ORCA-II ERG	
C4742-98-EWG(2)	ORCA-II EWG	
C4742-98-2xKAG(2) / -2xKWG(2)	ORCA-II BTG 1024	
C4742-98-2xLAG(2) / -2xLWG(2)	ORCA-II BTG 512	
C8484-xxG		(*2)
C8484-xxG01		
C9440-xxG		(*2)

Active Silicon Phoenix

DCAM Version		
DCAM Module	3.00.352.3032 2.1.3.3032	(for C9100 and C7780) (for ORCA, ORCA II, C7190)
DRIVER	5.32.0.3032	
Supported Board		
AS-PHX-D24CL-PCI32	Phoenix Std CameraLink	(for PowerMac)
AS-PHX-D24CL-PE1	Phoenix PCIe CameraLink	(for IntelMac)
AS-PHX-D36-PCI64U	Phoenix Pro LVDS	(for PowerMac)
AS-PHX-D36-PE1	Phoenix PCIe LVDS	(for IntelMac)
AS-PHX-D48CL-PCI64U	Phoenix Pro CameraLink	(for PowerMac)
AS-PHX-D48CL-PE1	Phoenix PCIe CameraLink	(for IntelMac)
Supported Camera		
C4742-95-12NR(B)	ORCA-NR	
C4742-95-12ER	ORCA-ER	
C4742-98-2xER	ORCA-II ER	
C4742-98-2xKA(2) / -2xKW(2)	ORCA-II BT 1024	
C4742-98-2xLA(2) / -2xLW(2)	ORCA-II BT 512	
C7190-5x		(*3)
C7780-xx		
C9100-01/02/11/12		
C9100-13	ImagEM	
C9100-14	ImagEM 1K	

Notes:

*New: New supported camera.

*1: CRi LCD filter driving is supported.

*2: These camera should not be connected to Mac's on-board Firewire port because it supplies about 30V which may damage the camera. Please use a PCI Firewire Adapter card.

*3: Special Digital I/F cable for Phoenix is required:

(a) HPK cable part number: A9608-95 (Digital I/F cable for Phoenix/C7190-20)

or

(b) For C7190-2x/4x/5x, HC/ASL cable part number:

WB3021 (CABLE FOR DIG-36-PE1/DIG-36 PHOENIX BOARD AND C7190-2x/4x/5x, ORCA-100/II/ER)

or

(c) For C4742-95-12HR, HC/ASL cable part number:

WB3021-HR (CABLE FOR DIG-36-PE1/DIG-36 PHOENIX BOARD AND ORCA-HR)

with Phoenix Serial Port jumpers set for RS232 for MacOSX.

All LVDS/RS-422 cameras works with Active Silicon Phoenix LVDS

All CamLink cameras work with Active Silicon Phoenix CamLink